


Mess.-Nr. 58923 **Datum** Donnerstag, 13 Dezember 2012 03:18 **Bearbeitungs- Nr.:**
Material Protein **S120218**
Bemerkung wetted with 4-5drops of tensid-solution (= 1drop of tensid: Dusazin 901 in 10ml of water), + 10ml of water and 30s ultrasonic bath LABORETTE 17

Berechnung Fraunhofer broad **TradeOff** 1,000000e+003
Scans 100 **Scans** 100 **Messkanäle** 102 **Strahlabsorption** 13,00 %
Feinbereich **Grobbereich**
Messbereich 0,08 [μm] - 2000,00 [μm] **Pumpe** 5 **Ultraschall** 10

%	< μm
5	0,78
10	1,22
15	1,74
20	2,39
25	3,28
30	4,52
40	7,92
50	11,26
60	14,44
70	17,88
80	22,04
90	28,17
95	33,41
98	39,63
99	43,74

%	< μm
0,0	0,1
1,9	0,5
7,6	1,0
17,1	2,0
31,6	5,0
46,1	10,0
75,4	20,0
92,1	30,0
98,1	40,0
99,7	50,0
100,0	60,0
100,0	70,0
100,0	80,0
100,0	90,0
100,0	100,0

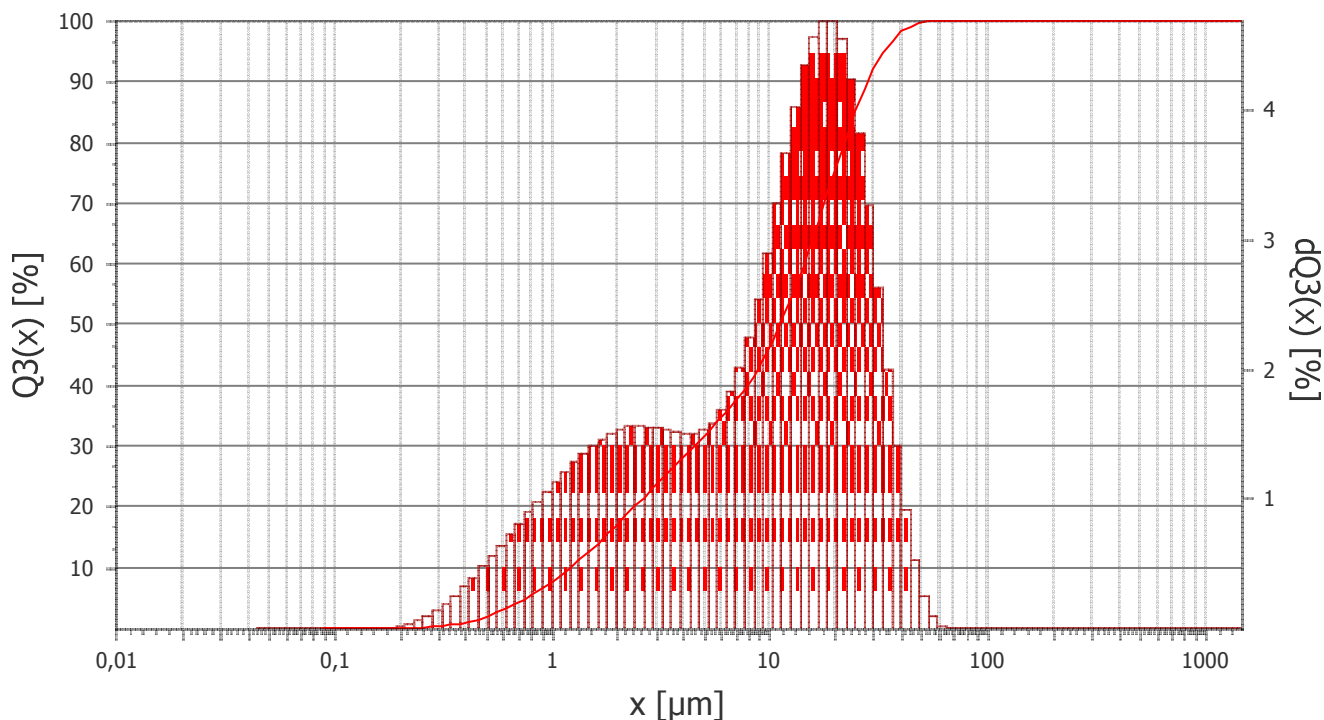

 58923 dQ3(x) — 58923 Q3(x)